

NLDFT Model Downloading

In addition to the full list of Micromeritics NLDFT models, a link to download new models is provided so that Micromeritics users may easily upgrade their software to include the latest innovations in isotherm modeling.

Instructions for downloading models

1. After logging into your customer portal, choose the [Download](#) link located in the right column.
 2. Click on the Model of your choice.
- Have the new model in the appropriate Micromeritics models directory as indicated below.

Micromeritics Application	Default model directory
AWM DataMaster	C:\DATAMSTR\MODELS
ASAP 2020	C:\2020\MODELS
ASAP 2420	C:\2420\MODELS
ASAP 3020	C:\WIN3020\MODELS
ASAP 2390	C:\WIN2390\MODELS

Restart your application and the new models are available via the DFT reports.